


Moisture analyser KERN MLS-A



Moisture analyser with 49 memories for individual drying processes

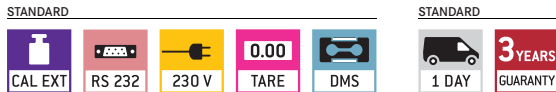

KERN	MLS 65-3A	MLS 150-2A
Readout [d]	0,001 g / 0,01 %	0,01 g / 0,05 %
Weighing range [Max]	65 g	150 g
Reproducibility, weight of sample 2 g	0,5 %	-
Reproducibility, weight of sample 10 g	0,05 %	0,2 %
Display after drying (Display can be switched over at any time)		
Moisture [%] = Weight loss (WL) from SW (startweight)	0 - 100 %	
Dry mass [%] = Residual weight (RW) from SW	100 - 0 %	
ATRO [%] [(SW - RW) : RW] · 100%	0 - 999 %	
Residual weight [g]	Absolute value in [g]	
Temperature range	50° - 160°C in steps up to 1°C	
Drying modes	<ul style="list-style-type: none"> • Standard drying • Drying in levels • Gentle drying • Pre-heat level, can be switched on 	
Switch off criteria	<ul style="list-style-type: none"> • When the set time has expired 3 min - 99 min • When the weight loss per unit of time falls below the target value (both values freely adjustable mg/sec) 	
Recall of measurement	Interval adjustable from 1 sec - 99 sec (Also in combination with KERN YKB-01N printer or PC)	
Overall dimensions WxDxH	250x360x175 mm	
Net weight	approx. 6 kg	

New model
























Features

- LCD display, digit height 25 mm
- 1 Previous drying time
- 2 Current temperature
- 3 Moisture display in %
- 4 Current moisture content in %
- 5 Halogenquartz glass heater 400 W
- The last value measured remains on the display until it is replaced by a new measurement
- 6 10 sample plates included
- Table of applications: There are many practical examples in the operating instructions
- Memory for automatic sequence of 49 complete drying processes. 10 free digits for each memory
- Easy to use
- Display in German and English

Options

- 6 Sample plates aluminium, Ø 92 mm. Unit of 80 pieces, **KERN MLB-A01**
- Round fiberglass filter e.g. for samples that splash or become encrusted. Unit of 80 pieces, **KERN RH-A02**
- Temperature calibration set consists of measuring sensor and display device, **KERN MLS-A04**
- Standard printer, **KERN YKB-01N**

KERN Pictograms

 Internal adjusting (CAL): Quick setting of the balance's accuracy with internal adjusting weight (motordriven).	 Adjusting program (CAL): For quick setting of the balance's accuracy. External adjusting weight required.	 Weighing units: Can be switched to e.g. non-metric units at the touch of a key. See balance model. Please refer to KERN's website for more details.
 Rechargeable battery pack: rechargeable set.	 Power supply: integrated in balance. 230V/50Hz in Germany. On request GB or USA version.	 Mains adapter: 230V/50Hz in standard version for Germany. On request GB or USA version.
 Battery operation: The battery type is specified for each device.	 Piece counting: Reference quantities selectable. Display can be switched from piece to weight	 Spray and dust protection IPxx The type of protection is shown by the pictogram. For details see the glossary.
 Data interface: The type of interface is shown by the pictogram. See the glossary for further details.	 GLP/ISO record keeping of weighing data with date, time and identification-no. Only with printers from KERN.	 Tare: put back displayed value to "0" during weighing process, e.g. to add or remove weighing material from a container.
 Animal weighing program: Strong vibrations are filtered out to determine a stable weighing result.	 Stainless steel: the balance is protected against corrosion.	 Warranty: The warranty period is shown in the pictogram.
 Net-total weighing: weight of tare cup and weight of components memorized in two separate stores.	 Suspended weighing: load support with hook on the underside of the balance.	 Verification possible. The time required for verification is specified in the pictogram.
 Percentage determination: Displays the deviation from the reference weight (100%) in % instead of grams.	 Weighing with tolerance range: Upper and lower limiting can be programmed individually, e.g. dosing/sorting.	 DKD calibration possible. The time required for DKD calibration is specified in the pictogram.
 Package shipment via courier.	 Pallet shipment via freight forwarder.	[DMS] For details on weighing technologies, see the glossary.

Precision is our business

To ensure the high level of precision of your balance, KERN offers the appropriate test weight package for your balance. This consists of the test weight, box and DKD calibration certificate, as proof of its accuracy. The best way to ensure proper balance calibration.

In the extensive KERN test weight range, you will find test weights in the international OIML error limit classes: E1, E2, F1, F2, M1, M2, and M3 with weights from 1 mg to 2000 kg.

The KERN DKD calibration laboratory for electronic balances and weights has been accredited by DKD since 1994 and today is one of the most modern and best-equipped DKD calibration laboratories for balances, test weights and force-measurement in Europe.
(DKD = German Calibration Service)

Thanks to the high level of automation, we can carry out DKD calibration of balances, test weights and force-measuring devices 24 hours a day, 7 days a week.

Range of services:

- DKD calibration of balances with a maximum load of up to 6 t
- DKD calibration of weights in the range of 1 mg – 500 kg
- Database supported management of checking equipment and reminder service
- Calibration of force-measuring devices
- DKD calibration certificates in the following languages
D, GB, F, I, E, NL, PL

Do you have questions about your scale, the corresponding test weight or the calibration service? Your KERN specialist dealer will be pleased to assist you.

Your KERN specialist dealer:

KERN – Professional measuring. Measuring technology and testing services from a single source



German Excellence Group
Member